

## CleanMax® Clean Manufactured Sterile Apparel – Donning Procedure

- Good Hygiene should be maintained prior to coming to work.
- Remove all jewelry and personal items.



Don bouffant cap, be sure all hair is contained and ears are covered.



Don disposable shoe covers making sure entire surface of street shoe is contained.



Wash or sanitize hands. Make sure they dry completely.



Don first pair of sterile gloves without touching outside surface area of glove.



Sanitize gloves – this should be done after each step in the gowning process.



Don Hood – use ties to ensure proper fit.



Don Mask – adjust for proper fit (no gaps)



Don Coverall with approved aseptic technique. Be careful removing coverall from package. Touch only the inside of the coverall, do not let coverall touch operator, floor or any surface.



Don first boot cover with approved aseptic technique, then place first donned boot onto "clean side" of gowning area. Repeat with second boot cover.



Don Goggles – adjust for sealed fit.



Don Second Pair of Sterile Gloves, completely covering the first pair.



Do final check in mirror prior to entering sterile cleanroom.

## CleanMax® Clean Manufactured Sterile Apparel – Doffing Procedure

- Exit sterile cleanroom. Remove items in designated area in reverse order of gowning and discard.



Remove outer gloves.



Remove goggles.



Remove shoe covers.



Remove mask.



Remove coverall.



Remove hood.



Remove inner gloves.



Remove shoe covers.



Remove bouffant cap.



800-645-9291 | 256-350-3873 | [info@lakeland.com](mailto:info@lakeland.com) | [lakeland.com/cleanmax](http://lakeland.com/cleanmax)

Warning: Cleanroom apparel should not be used around heat, flames, sparks or in potentially flammable or explosive environments.  
Cleanroom fabrics should have slip-resistant materials on the outer sole of boots, shoe covers, or other garment surfaces in conditions where slipping could occur.  
4311/6-17-2020/